


NIRLAB

Quick Start Manual

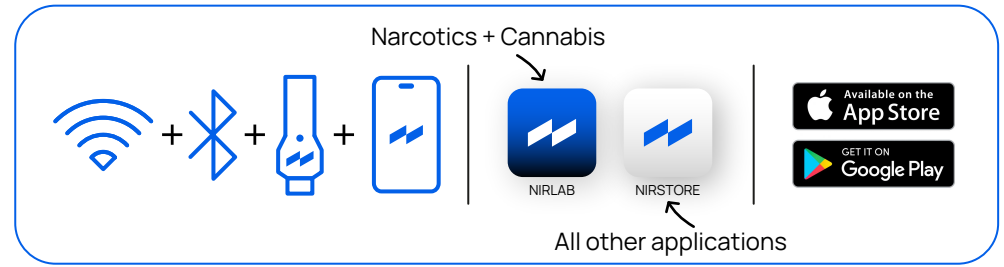

Get started

1. Login on the APP
2. Click on the device button
3. Pair

Calibrate

1. Click the calibration button if orange
2. Calibration type R99-mirror Calibration accessory NO ACCESSORY
3. Place the device on the calibration mirror Calibration done!

Scan tips

Direct contact

Point downwards

Aluminium cup

Clean w/ ethanol

Sample Scan

Sample information / Purity average
Multiple scans for 1 sample

1. Click the sample scan button
2. Name Attributes Photo Sample info
3. Your Substance % Purity average

Generate report (PDF)

Quick Scan

Generates automatic name
1 Analysis only

1. Click the quick scan button
2. Your Substance Purity %

Generate report (PDF)

Spectrum

or click only!

Turn off & more

1. Click for 3 Seconds Or

Narcotics & Cannabis : <https://app.nirlab.com>

See more here

All other applications: <https://nirstore.nirlab.com>